


PRISM MICROSCOPE

Acorn Naturalists #T-15293

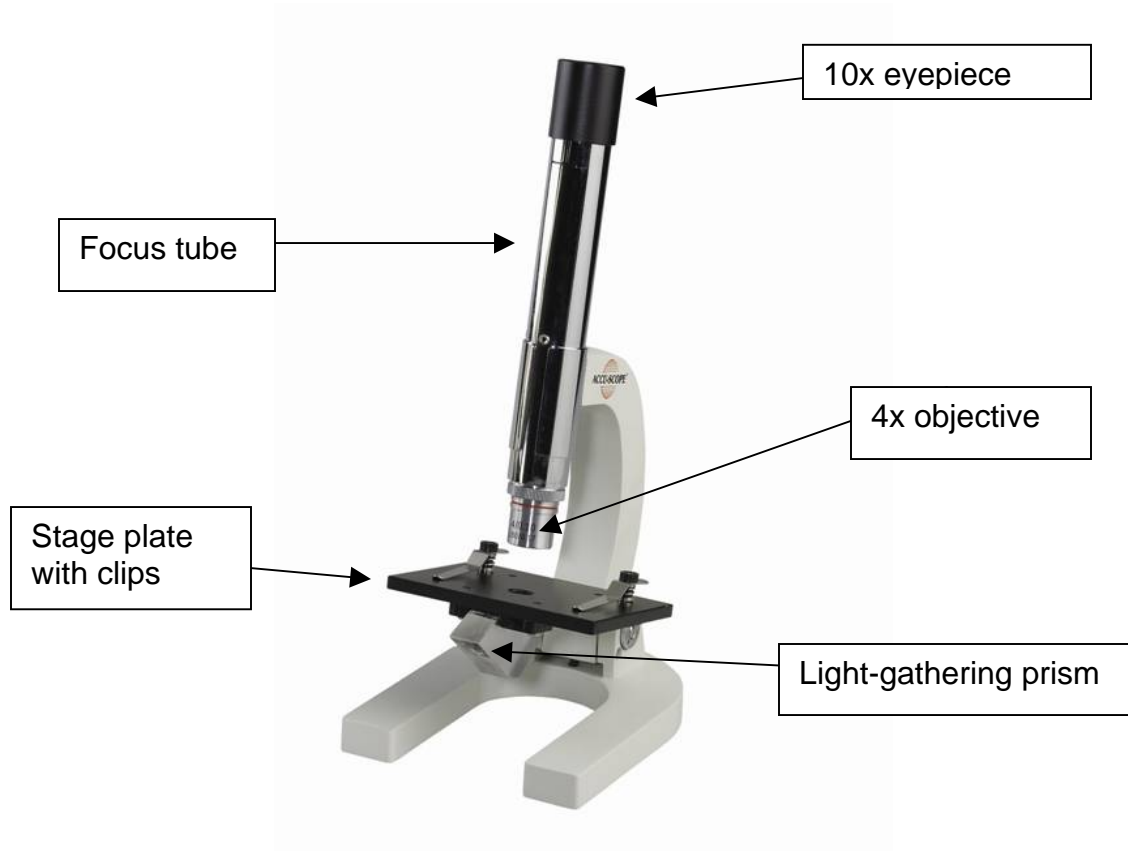

Prism microscope information:

- This instrument has a 4x objective and uses ambient or room light.
- The working distance of the 4x objective is approximately 1.6mm.
- To view a slide or other item, place on the stage plate, securing with the clips and simply slide the focus tube up or down until the object comes into focus.
- When needing a finer adjustment, rotate the focus tube clockwise or counter clockwise.
- Lenses may be cleaned occasionally with lens paper and mild lens cleaner such as used for eyeglasses. No other maintenance is needed.